

# Product information



Information about other products is available at: [www.IBL-America.com](http://www.IBL-America.com)

## ImmunoGuide®

### Instructions for Use

## Quality Control (QC) Sera

For Adalimumab ELISA coded TM09008

**REF: TM09008-QC**



2 Vials



2-8°C

**For Research Use Only - Not for Use in Diagnostic Procedures**

Manufactured for  
**Immuno-Biological Laboratories Inc. (IBL-America)**  
8201 Central Avenue, NE, Suite P  
Minneapolis, MN 55432  
Tel: 763-780-2955 Toll Free: 1-888-523-1246  
[www.IBL-America.com](http://www.IBL-America.com)

## 1. Product Description

**Catalogue Number:** TM09008-QC

**Size:** 2 vials, lyophilized

**Adalimumab-High Control Lyophilized (1 vial containing 0.5mL):** Human serum containing the drug Adalimumab at 5 µg/mL with added stabilizer and preservative.

**Adalimumab-Low Control Lyophilized (1 vial containing 0.5mL):** Human serum containing the drug Adalimumab at 1 µg/mL with added stabilizer and preservative.

**Application:**Used as a complementary component for TM09008 & TM09009 ELISAs.

**Safety Precautions:** The human sera contained in this product have been tested and found negative for HCV, HBsAg and HIV 1&2 by ELISA methods. Since no one test is 100% accurate, the product should be handled as if potentially infectious. Operators should wear gloves and protective clothing when handling these products. These products contains 0.09% sodium azide as preservative and is classified as non-hazardous according to EC directives.

## 2. STORAGE AND RECONSTITUTION INSTRUCTIONS

This lyophilized Quality Control (QC) material is shipped at ambient temperature. Upon receipt the lyophilized product should be stored at 2-8 °C and the non-reconstituted vial can be used until the expiry date as indicated on the label. After reconstitution the product should be stored at 2-8 °C and has a shelf life of 1 week. Storing of reconstituted products at  $\leq -50^{\circ}\text{C}$  prolongs their shelf life up to 8 months.

**2.1.** Reconstitute each vial, containing lyophilized powder with 0.5 mL of distilled water.

**2.2.** After addition of distilled water, recap the vial and allow the reconstituted serum to equilibrate for at least 30 minutes before mixing.

**2.3.** Mix the reconstituted serum by cautiously turning up and down 15-20 times. Do not vortex or mix vigorously.