

Product information



Information about other products is available at: www.IBL-America.com

ImmunoGuide®

Instructions for Use

Quality Control (QC) Sera

For Infliximab ELISA coded TM09001 and TM09002

REF: TM09001-QC



2 Vials



2-8°C

For Research Use Only - Not for Use in Diagnostic Procedures

Manufactured for
Immuno-Biological Laboratories Inc. (IBL-America)
8201 Central Avenue, NE, Suite P
Minneapolis, MN 55432
Tel: 763-780-2955 Toll Free: 1-888-523-1246
www.IBL-America.com

1. Product Description

Catalogue Number: TM09001-QC

Size: 2 vials, lyophilized

Infliximab-High Control Lyophilized: Human serum containing the drug Infliximab at 10 µg/mL with added stabilizer and preservative (1 vial containing 0.5 mL).

Infliximab-Low Control Lyophilized: Human serum containing the drug Infliximab at 1 µg/mL with added stabilizer and preservative (1 vial containing 0.5 mL).

Application: Used as a complementary component for TM09001 and TM09002 ELISAs.

Safety Precautions: The human sera contained in this product have been tested and found negative for HCV, HBsAg and HIV 1&2 by ELISA methods. Since no one test is 100% accurate, the product should be handled as if potentially infectious. Operators should wear gloves and protective clothing when handling these products.

2. STORAGE AND RECONSTITUTION INSTRUCTIONS

This lyophilized Quality Control (QC) material is shipped at ambient temperature. Upon receipt the lyophilized product should be stored at 2-8 °C and the non-reconstituted vial can be used until the expiry date as indicated on the label. After reconstitution the product should be stored at 2-8 °C and has a shelf life of 1 week. Storing of reconstituted products at ≤ -50°C prolongs their shelf life up to 8 months.

2.1. Reconstitute each vial, containing lyophilized powder with 0.5 mL of distilled water.

2.2. After addition of distilled water, recap the vial and allow the reconstituted serum to equilibrate for at least 30 minutes before mixing.

2.3. Mix the reconstituted serum by cautiously turning up and down 15-20 times. Do not vortex or mix vigorously.