

ORGENTEC Diagnostika GmbH

Carl-Zeiss-Straße 49-51

55129 Mainz - Germany

Phone: +49 (0) 61 31 / 92 58-0

Fax: +49 (0) 61 31 / 92 58-58

Internet: www.orgentec.com



**For research use only.
Not for diagnostic procedures.**



RUO 601 Anti-CCP hs® (high sensitive)

INTENDED USE

Anti-CCP hs® (high sensitive) is an ELISA test system for the quantitative measurement of IgG class autoantibodies against cyclic citrullinated peptides (CCP) in human serum or plasma. This product is intended for research use only. Not for diagnostic procedures.



Distributed By:

IBL-America, Inc.

8201 Central Ave NE, Suite P

Minneapolis, MN 55432, USA

info@ibl-america.com

(888) 523 1246

SYMBOLS USED ON LABELS



Manufacturer



Catalogue number



Sufficient for 96 determinations



Batch code



Use by



Temperature limitation



Keep away from sunlight



Do not reuse



Date of manufacture



Consult instructions for use

MICROPLATE

Microplate

CALIBRATOR A

Calibrator

CALIBRATOR B

Calibrator

CALIBRATOR C

Calibrator

CALIBRATOR D

Calibrator

CALIBRATOR E

Calibrator

CALIBRATOR F

Calibrator

CONTROL +

Control positive

CONTROL -

Control negative

DILUENT

Sample Buffer P

CONJUGATE

Enzyme Conjugate

TMB

TMB Substrate

STOP

Stop solution

WASH

Wash Buffer

RTU

Ready to use

PRINCIPLE OF THE TEST

Highly purified cyclic citrullinated vimentin peptides (CCP) is bound to microwells.

The determination is based on an indirect enzyme linked immune reaction with the following steps:

Specific antibodies in the sample bind to the antigen coated on the surface of the reaction wells. After incubation, a washing step removes unbound and unspecifically bound serum or plasma components. Subsequently added enzyme conjugate binds to the immobilized antibody-antigen-complexes. After incubation, a second washing step removes unbound enzyme conjugate. After addition of substrate solution the bound enzyme conjugate hydrolyses the substrate forming a blue coloured product. Addition of an acid stops the reaction generating a yellow end-product. The intensity of the yellow color

correlates with the concentration of the antibody-antigen-complex and can be measured photometrically at 450 nm.

WARNINGS AND PRECAUTIONS

- All reagents of this kit are intended for professional use only.
- Components containing human serum were tested and found negative for HBsAg, HCV, HIV1 and HIV2 by FDA approved methods. No test can guarantee the absence of HBsAg, HCV, HIV1 or HIV2, and so all human serum based reagents in this kit must be handled as though capable of transmitting infection.
- Bovine serum albumin (BSA) used in components has been tested for BSE and found negative.
- Avoid contact with the substrate TMB (3,3',5,5'-Tetramethyl-benzidine).
- Stop solution contains acid, classification is non-hazardous. Avoid contact with skin.
- Control, sample buffer and wash buffer contain sodium azide 0.09% as preservative. This concentration is classified as non-hazardous.
- Enzyme conjugate contains ProClin 300 0.05% as preservative. This concentration is classified as non-hazardous.

During handling of all reagents, controls and serum samples observe the existing regulations for laboratory safety regulations and good laboratory practice:








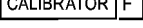
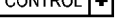


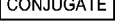
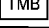
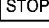


- First aid measures: In case of skin contact, immediately wash thoroughly with water and soap. Remove contaminated clothing and shoes and wash before reuse. If system fluid comes into contact with skin, wash thoroughly with water. After contact with the eyes carefully rinse the opened eye with running water for at least 10 minutes. Get medical attention if necessary.
- Personal precautions, protective equipment and emergency procedures:

Observe laboratory safety regulations. Avoid contact with skin and eyes. Do not swallow. Do not pipette by mouth. Do not eat, drink, smoke or apply makeup in areas where specimens or kit reagents are handled. When spilled, absorb with an inert material and put the spilled material in an appropriate waste disposal.

- Exposure controls / personal protection: Wear protective gloves of nitril rubber or natural latex. Wear protective glasses. Used according to intended use no dangerous reactions known.
- Conditions to avoid: Since substrate solution is light-sensitive. Store in the dark.
- For disposal of laboratory waste the national or regional legislation has to be observed.

Observe the guidelines for performing quality control in medical laboratories by assaying control sera.

CONTENTS OF THE KIT

RUO 601	 96	Sufficient for 96 determinations
	1	One divisible microplate consisting of 12 modules of 8 wells each. Ready to use. Product code on module: CCP
 A	1.5 ml	Calibrator A 0 U/ml, containing serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use.
 B	1.5 ml	Calibrator B 20 U/ml, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use.
 C	1.5 ml	Calibrator C 40 U/ml, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use.
 D	1.5 ml	Calibrator D 100 U/ml, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use.
 E	1.5 ml	Calibrator E 300 U/ml, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, NaN3 0.09%), yellow. Ready to use.
 F	1.5 ml	Calibrator F 1000 U/ml, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use.
 +	1.5 ml	Control positive, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use. The concentration is specified on the certificate of analysis.
 -	1.5 ml	Control negative, containing CCP antibodies in a serum/buffer matrix (PBS, BSA, detergent, NaN3 0.09%), yellow. Ready to use. The concentration is specified on the certificate of analysis.
	20 ml	Sample Buffer P, containing PBS, BSA, detergent, preservative sodium azide 0.09%, yellow, concentrate (5 x).
	15 ml	Enzyme Conjugate containing anti-human IgG antibodies, HRP labelled; PBS, BSA, detergent, preservative PROCLIN 0.05%, light red. Ready to use.
	15 ml	TMB Substrate; containing 3,3', 5,5'- Tetramethylbenzidin, colorless. Ready to use.
	15 ml	Stop solution; contains acid. Ready to use.
	20 ml	Wash Buffer, containing Tris, detergent, preservative sodium azide 0.09%; 50 x conc.
	1	Certificate of Analysis

MATERIALS REQUIRED

- Microplate reader capable of endpoint measurements at 450 nm; optional: reference filter at 620 nm
- Data reduction software
- Multi-channel dispenser or repeatable pipette for 100 µl
- Vortex mixer
- Pipettes for 10 µl, 100 µl and 1000 µl
- Laboratory timing device
- Distilled or deionised water
- Measuring cylinder for 1000 ml and 100 ml
- Plastic container for storage of the wash solution

This ELISA assay is suitable for use on open automated ELISA processors. Each assay has to be validated on the respective automated system. Detailed information is provided upon request.

SPECIMEN COLLECTION, STORAGE AND HANDLING

- Collect whole blood specimens using acceptable medical techniques to avoid hemolysis.
- Allow blood to clot and separate the serum or plasma by centrifugation.
- Test serum should be clear and non-hemolyzed. Contamination by hemolysis or lipemia should be avoided, but does not interfere with this assay.
- Specimens may be refrigerated at 2-8°C for up to five days or stored at -20°C up to six months.
- Avoid repetitive freezing and thawing of serum or plasma samples. This may result in variable loss of antibody activity.
- Testing of heat-inactivated sera is not recommended.

STORAGE AND STABILITY

- Store test kit at 2-8°C in the dark.
- Do not expose reagents to heat, sun, or strong light during storage and usage.
- Store microplate sealed and dessicated in the clip bag provided.
- Shelf life of the unopened test kit is 18 months from day of production.
Unopened reagents are stable until expiration of the kit. See labels for individual batch.
- Diluted Wash Buffer and Sample Buffer are stable for at least 30 days when stored at 2-8°C.
We recommend consumption on the same day.

PROCEDURAL NOTES

- Do not use kit components beyond their expiration dates.
- Do not interchange kit components from different lots and products.
- All materials must be at room temperature (20-28°C) prior to use.
- Prepare all reagents and samples. Once started, perform the test without interruption.
- Double determinations may be done. By this means pipetting errors may become obvious.
- Perform the assay steps only in the order indicated.
- Always use fresh sample dilutions.
- Pipette all reagents and samples into the bottom of the wells.
- To avoid carryover or contamination, change the pipette tip between samples and different kit controls.
- Wash microwells thoroughly and remove the last droplets of wash buffer.
- All incubation steps must be accurately timed.
- Do not re-use microplate wells.

PREPARATION OF REAGENTS

WASH

Dilute the contents of one vial of the buffered wash solution concentrate (50x) with distilled or deionised water to a final volume of 1000 ml prior to use.

DILUENT

Sample Buffer P: Prior to use dilute the contents (20 ml) of one vial of sample buffer 5x concentrate with distilled or deionised water to a final volume of 100 ml.

Preparation of samples

Dilute samples 1:100 before the assay: Put 990 µl of prediluted sample buffer in a polystyrene tube and add 10 µl of sample. Mix well. Note: Calibrators / Controls are ready to use and need not be diluted.

TEST PROCEDURE

Prepare enough microplate modules for all calibrators / controls and samples.

1. Pipette 100 µl of calibrators, controls and prediluted samples into the wells.
Incubate for 30 minutes at room temperature (20-28 °C).
Discard the contents of the microwells and wash 3 times with 300 µl of wash solution.
2. Dispense 100 µl of enzyme conjugate into each well.
Incubate for 15 minutes at room temperature.
Discard the contents of the microwells and wash 3 times with 300 µl of wash solution.
3. Dispense 100 µl of TMB substrate solution into each well.
Incubate for 15 minutes at room temperature
4. Add 100 µl of stop solution to each well of the modules
Incubate for 5 minutes at room temperature.
Read the optical density at 450 nm (reference 600-690nm) and calculate the results.
The developed colour is stable for at least 30 minutes. Read during this time.

Example for a pipetting scheme:

	1	2	3	4	5	6	7	8	9	10	11	12
A	A	P1										
B	B	P2										
C	C	P3										
D	D											
E	E											
F	F											
G	C+											
H	C-											

P1, ... sample A-F calibrators C+, C- controls

VALIDATION

Test results are valid if the optical densities at 450 nm for calibrators / controls and the results for controls comply with the reference ranges indicated on the Certificate of Analysis enclosed in each test kit.
If these quality control criteria are not met the assay run is invalid and should be repeated.

CALCULATION OF RESULTS

For quantitative results plot the optical density of each calibrator versus the calibrator concentration to create a calibration curve. The concentration of samples may then be estimated from the calibration curve by interpolation.
Using data reduction software a 4-Parameter-Fit with lin-log coordinates for optical density and concentration is the data reduction method of choice.

ASSAY CHARACTERISTICS

Calibration

This assay system is calibrated in relative arbitrary units. It is calibrated against an external anti-CCP Assay, since no international reference sera for RA diagnostic are available so far.

Measuring range

The calculation range of this ELISA assay is 0 - 1000 U/ml

Interfering substances

No interference has been observed with haemolytic (up to 1000 mg/dl) or lipemic (up to 3 g/dl triglycerides) sera or plasma, or bilirubin (up to 40 mg/dl) containing sera or plasma. Nor have any interfering effects been observed with the use of anticoagulants (Citrate, EDTA, Heparine). However for practical reasons it is recommended that grossly hemolyzed or lipemic samples should be avoided.

LIMITATIONS OF THE PROCEDURE

For research use only. Not for diagnostic procedures.

REFERENCES

1. Bang H, Egerer K, Gauliard A, et al. Mutation and citrullination modifies vimentin to a novel autoantigen for rheumatoid arthritis. *Arthritis Rheum.* 56(8):2503, 2007.
2. Aletaha D, et al. Rheumatoid arthritis classification criteria: An American College of Rheumatology/European League Against Rheumatism collaborative initiative. *Arthritis & Rheumatism* 62 (9):2569-2581, 2010.
3. Pruijn GJ, Wiik A, van Venrooij WJ. The use of citrullinated peptides and proteins for the diagnosis of rheumatoid arthritis. *Arthritis Res Ther* 12 (1):203, 2010.
4. Snir O, Widhe M, Hermansson M, von Spee C, Lindberg J, Hensen S, Lundberg K, Engstrom A, Venables PJ, Toes RE, Holmdahl R, Klareskog L, Malmstrom V. Antibodies to several citrullinated antigens are enriched in the joints of rheumatoid arthritis patients. *Arthritis Rheum* 62 (1):44-52, 2009.
5. Van Steendam K, Tilleman K, Deforce D. The relevance of citrullinated vimentin in the production of antibodies against citrullinated proteins and the pathogenesis of rheumatoid arthritis. *Rheumatology (Oxford)*, 2011.
6. Van Steendam K, Tilleman K, De Ceuleneer M, De Keyser F, Elewaut D, Deforce D. Citrullinated vimentin as an important antigen in immune complexes from synovial fluid of rheumatoid arthritis patients with antibodies against citrullinated proteins. *Arthritis Res Ther* 12 (4):R132, 2010.
7. Syversen SW, Goll GL, van der Heijde D, Landewe R, Lie BA, Odegard S, Uhlig T, Gaarder PI, Kvien TK. Prediction of radiographic progression in rheumatoid arthritis and the role of antibodies against mutated citrullinated vimentin: results from a ten-year prospective study. *Ann.Rheum.Dis.* 69 (2):345-351, 2009.
8. Innala L, Kokkonen H, Eriksson C, Jidell E, Berglin E, Dahlqvist SR. Antibodies against mutated citrullinated vimentin are a better predictor of disease activity at 24 months in early rheumatoid arthritis than antibodies against cyclic citrullinated peptides. *J Rheumatol.* 35(6):1002, 2008.
9. Mathsson L, Mullazehi M, Wick MC, et al. Antibodies against citrullinated vimentin in rheumatoid arthritis: higher sensitivity and extended prognostic value concerning future radiographic progression as compared with antibodies against cyclic citrullinated peptides. *Arthritis Rheum.* 58(1):36, 2008.
10. Soós L, Szekanecz Z, Szabó Z, et al. Clinical evaluation of anti-mutated citrullinated vimentin by ELISA in rheumatoid arthritis. *J Rheumatol.* 34(8):1658, 2007.
11. Bizzaro N, Tonutti E, Tozzoli R, Villalta D. Analytical and diagnostic characteristics of 11 2nd- and 3rd-generation immunoenzymatic methods for the detection of antibodies to citrullinated proteins. *Clin Chem.* 53 (8):1527, 2007.
12. Coenen D, Verschueren P, Westhovens R, Bossuyt X. Technical and diagnostic performance of 6 assays for the measurement of citrullinated protein/peptide antibodies in the diagnosis of rheumatoid arthritis. *Clin Chem.* 53 (3):498, 2007.
13. Dejaco C, Klotz W, Larcher H, Duftner C, Schirmer M, Herold M. Diagnostic value of antibodies against a modified citrullinated vimentin in rheumatoid arthritis. *Arthritis Res Ther.* 8(4):R119, 2006.
14. Keskin G, Inal A, Keskin D, et al. Diagnostic utility of anti-cyclic citrullinated peptide and anti-modified citrullinated vimentin antibodies in rheumatoid arthritis. *Protein Pept Lett.* 15(3):314 2008.
15. Poulosom H, Charles PJ. Antibodies to citrullinated vimentin are a specific and sensitive marker for the diagnosis of rheumatoid arthritis. *Clin Rev Allergy Immunol.* 34(1):4 2008.
16. Song JS, Park GB, Park AJ. Comparison of anti-mutated citrullinated vimentin with anti-cyclic citrullinated peptide and rheumatoid factors for the diagnostic value of rheumatoid arthritis [in Korean]. *J Korean Rheum Assoc.* 14(3):235 2007.
17. Ursum J, Nielen MM, van Schaardenburg D, et al. Antibodies to mutated citrullinated vimentin and disease activity score in early arthritis: a cohort study. *Arthritis Res Ther.* 10(1):R12, 2008.
18. Besada E, Nikolaisen C, Nossent H. Diagnostic value of antibodies against mutated citrullinated vimentin for rheumatoid arthritis. *Clin Exp Rheumatol.* 29(1):85 2011.

Change Control

Former version: -

Reason for revision: *first edition*